

SOFTWARE RELEASE NOTES

Experiment Manager v2.3.1

About this Software

Experiment Manager v2.3.1 is designed to run the LeviCell® 1.0 system. The Experiment Manager software is pre-installed on the Windows 10 Pro 64-bit computer provided with the LeviCell 1.0 instrument.

For use on PC only.

Item	Description
For Use With	LeviCell 1.0 System
Last Released Version	v2.2.1
New Software Version	v2.3.1
New Software Release Date	March 2025
New Features	<ul style="list-style-type: none">• Optimizes the image saving scheme and increases run storage capacity by 4X.
Bug Fixes and Improvements	<ul style="list-style-type: none">• Reduces the live fluorescence image frequency to prevent cell movement during levitation• Displays the split line during the levitation when it is pre-set• Limit the number of slow advances to 3 to prevent user error• General UI improvements

Experiment Manager Update Procedure

1. Go to www.levitasbio.com/support.

2. Download the **LeviCell Experiment Manager v2.3.1.1150.msi** file for the new software version to your LeviCell 1.0 computer.

If the computer is not connected to the internet, use an external storage device such as a USB drive to transfer the downloaded file to the LeviCell 1.0 computer.

3. Double-click to launch the .msi file.

4. A Windows protect screen will.

5. Confirm the publisher is LevitasBio, Inc., then select "**OK**" "**Run anyway**".

6. Follow the on-screen instructions to install the new software. *Upon completion, Windows may ask you to restart the computer.*

7. When installation is complete, a new shortcut for Experiment Manager v2.3 will appear on your desktop. To avoid confusion, be sure to remove shortcuts for earlier versions of Experiment Manager software.

8. To start Experiment Manager, double click on the software icon.



Contact Technical Support at support@levitasbio.com if any errors are encountered during the procedure.