

EOS Manager v3.0

About this Software

EOS Manager v3.0 is designed to run the LeviCell® EOS and LeviCell EOS TEC Systems. The EOS Manager software is pre-installed on the Windows 10 Pro 64-bit computer provided with LeviCell EOS Systems.

For use on LeviCell EOS Control PC only.

Item	Description
For Use With	LeviCell EOS and LeviCell EOS TEC Systems
Last Software Version	v2.1.3
New Software Version	v3.0.3
Software Release Date	December 2025
New Features	- Run storage capacity is increased by 10X with advanced image compression technique. - Add pagination to the run history view.

EOS Manager Update Procedure

1. Go to support website www.levitasbio.com/support
2. Choose LeviCell EOS product on the left.
3. Download the new software version *.msi file to the LeviCell EOS computer. If the LeviCell EOS computer is not connected to the internet, use an external storage device (such as a USB) to transfer the downloaded file to the computer.
4. Double-click on the “EOS Manager 3.0.3.1112.msi” file.
5. A Windows protection screen will appear. Click “More Info” in the text.
6. Confirm that the publisher is LevitasBio, Inc, and then click “Run Anyway” button.
7. Follow the instruction to install the new software until you have completed the installation.

Contact support@levitasbio.com if any errors are encountered during the procedure.