

LeviMetrics v3.1

About this Software

LeviMetrics™ v3.1 is designed to visualize and perform analysis on experiment runs from the LeviCell® EOS System and LeviCell® EOS TEC System. The LeviMetrics software can be installed on any Windows 10 or 11 Pro 64-bit , 32 GB RAM PC computer or can be added to the provided Control PC that is included with LeviCell EOS system.

Item	Description
For Use With	LeviCell EOS and LeviCell EOS TEC Systems
Last Released Version	v2.1.3 / v3.0.1 (Elite)
New Software Version	v3.1.1
New Software Release Date	December 2025
New Features	• Profile Analysis Enhancement (Elite only) Automatic spectral crosstalk correction. Visualize individual or all stain results. • Software licensing for seamless Elite upgrade. • Visualize and analyze runs with compressed images.

LeviMetrics Update Procedure

1. Go to support website www.levitasbio.com/support
2. Choose LeviMetrics product on the left.
3. Download the new software version *.msi file to the LeviCell EOS computer or a computer with Windows 10 or Windows 11. If the computer is not connected to the internet, use an external storage device (such as a USB) to transfer the downloaded file to the computer.
4. Double-click on the “**LeviMetrics 3.1.1.368.msi**” file.
5. A Windows protection screen may appear. Click “**More Info**” in the text.
6. Confirm that the publisher is LevitasBio, Inc, and then click “**Run Anyway**” button.
7. Follow the instruction to install the new software until you have completed the installation.

Contact support@levitasbio.com if any errors are encountered during the procedure.