

# LeviMetrics v2.0

## About this Software

LeviMetrics™ v2.0 is designed to visualize and perform analysis on experiment runs from the LeviCell® EOS system. The LeviMetrics software can be installed on any Windows 10 or 11 Pro 64-bit , 32 GB RAM PC computer or can be added to the provided Control PC that is included with LeviCell EOS system.

Item	Description
For Use With	LeviCell EOS System
Last Released Version	v1.1
New Software Version	v2.0
New Software Release Date	Jan 2024
New Features	<p><b>New User Interface</b></p> <p><b>Analysis Modes</b>            Run information            - Display run info entered by the user during the experiment            - Option to edit Levitation Agent Conc, Notes            - Tracking of updates (user and date/time)</p> <p>Run Review            Levitation Review            - View levitation of each sample            - View in fluorescence mode            - Screen capture or create movie</p> <p>Collection Review            - View collection of samples            - Screen capture or create movie</p> <p>Fractionation Analysis            - View fractionation % for all lanes            - View alternative split line values and fractionation % or separation factors</p> <p>Levitation Analysis            - View Relative Intensity Graph            - Simulate alternative Levitation Agent concentration</p>

## LeviMetrics Update Procedure

1. Go to support website [www.levitasbio.com/support](http://www.levitasbio.com/support)
2. Download the new software version \*.msi file to the LeviCell EOS computer or a computer with Windows 10 or Windows 11. If the computer is not connected to the internet, use an external storage device (such as a USB) to transfer the downloaded file to the computer.
3. Double-click on the “**Install LeviMetrics v2.0.msi**” file.
4. A Windows protection screen may appear. Click “**More Info**” in the text.
5. Confirm that the publisher is LevitasBio, Inc, and then click “**Run Anyway**” button.
6. Accept the LeviMetrics License Agreement by checking the “I accept the terms in the License Agreement” box and click Install.
7. Follow the instruction to install the new software until you have completed the installation.

**Windows may ask you to restart the computer after installation.**

Contact [support@levitasbio.com](mailto:support@levitasbio.com) or call Technical Support at +1 (650) 204-1185 if any errors are encountered during the procedure.