

EOS Manager v1.2

About this Software

EOS Manager v1.2 is designed to run the LeviCell® EOS System. The EOS Manager software is preinstalled on the Windows 10 Pro 64-bit computer provided with LeviCell EOS System.

For use on LeviCell EOS Control PC only.

Item	Description
For Use With	LeviCell EOS System
Last Software Version	v1.1
New Software Version	v1.2
Software Release Date	Jan 2024
New Features	<ul style="list-style-type: none">• Added Nuclei protocol• Updated Run Summary screen• Updated EOS-4 cartridge image

EOS Manager Update Procedure

1. Go to support website www.levitasbio.com/support
2. Download the new software version *.msi file to the LeviCell EOS computer. If the LeviCell EOS computer is not connected to the internet, use an external storage device (such as a USB) to transfer the downloaded file to the computer.
3. Double-click on the “Install EOS Manager v1.2.msi” file.
4. A Windows protection screen will appear. Click “More Info” in the text.
5. Confirm that the publisher is LevitasBio, Inc, and then click “Run Anyway” button.
6. Follow the instruction to install the new software until you have completed the installation.

Windows may ask you to restart the computer after installation.

Contact support@levitasbio.com if any errors are encountered during the procedure.