

## EOS Manager v2.1

### About this Software

EOS Manager v2.1 is designed to run the LeviCell® EOS and LeviCell EOS TEC Systems. The EOS Manager software is preinstalled on the Windows 10 Pro 64-bit computer provided with LeviCell EOS Systems.

*For use on LeviCell EOS Control PC only.*

Item	Description
For Use With	LeviCell EOS and LeviCell EOS TEC Systems
Last Software Version	v1.2
New Software Version	v2.1
Software Release Date	July 2024
New Features	<p><b>Compatibility with EOS TEC Core hardware</b></p> <ul style="list-style-type: none"> <li>• Updated nuclei protocol</li> <li>• Enable temperature control runs for nuclei protocol                             <ul style="list-style-type: none"> <li>• Cold, Cool, CRT (Controlled Room Temperature for cell runs)</li> </ul> </li> <li>• Added post-run protocol for drying system after a cold or cool run                             <ul style="list-style-type: none"> <li>• Added parallel collection option</li> </ul> </li> <li>• CRT temperature run for cell runs (Ambient is removed)</li> <li>• Added scheduled task Heat Dry at set time or prior to shutdown</li> </ul>

### EOS Manager Update Procedure

1. Go to support website [www.levitasbio.com/support](http://www.levitasbio.com/support)
2. Download the new software version \*.msi file to the LeviCell EOS computer. If the LeviCell EOS computer is not connected to the internet, use an external storage device (such as a USB) to transfer the downloaded file to the computer.
3. Double-click on the “**Install EOS Manager v2.1.msi**” file.
4. A Windows protection screen will appear. Click “**More Info**” in the text.
5. Confirm that the publisher is LevitasBio, Inc, and then click “**Run Anyway**” button.
6. Follow the instruction to install the new software until you have completed the installation.

**Windows may ask you to restart the computer after installation.**

**Contact [support@levitasbio.com](mailto:support@levitasbio.com) if any errors are encountered during the procedure.**